## Using a Microphone with Soundcard Oscilloscope

- 1. Sign in to the computer.
- 2. Turn on the microphone preamp.
- 3. When your desktop appears, plug in the microphone in to the computer's front headset jack.

Select "Mic In" and click on <OK>.







4. Start the Scope application.



5. If Scope does not react to the microphone:

Select the "Settings" tab and then

Input: Jack Mic (Realtek Audio).

6. If Scope still does not show a microphone signal, check the Windows Recording level.

Exit	Oscilloscope X-Y graph Frequency Signal generator Extras Settings
Amplitude CH1 Amplitude CH2 [1/Div] 10m 10m 10m 100m	Windows Sound Parameters           Open Audio Mixer         Audio Devices           Output         Output         Speakers / Headphones (Realtek million)           Image: Devices         Devices         Devices
tm 30m 0fiet 00m	auto ensible Data Format 44100 Sampleo/s 16 Bits/sample
Time (s) 100m 10m, 1	Scope Parameters On-screen values update period 500 ms Averaging time for f meaurement 500 ms
1m 10 Edge	Change Language Calibration of amplitude 10000 V/Dh Save and Restore Settings
Run/Stop	Save current settings Read settings from file Restore last settings Default file
Channel Mode	3

					- 6	X
	+ + + T 🛄 > Control Ranel > All	Control Parel Berry		D Seven Comm Pase )	Salect Central Panel	2
Control Panel   Sounds   Recording	Adjust your computer's settings			View by Lage core *	en by: Large store *	
Jack mic	Administrative lools	AutoPlay ® Sound	× 💐	BitLocker Drive Encryption	sker Drive Encryption	
<properties></properties>	Color Management	Configur Sated arcoard og	Search Communications denice bedgen to modelly do settings	Date and Time	and Time	
	Delault Programs	Dell Aud	t Delay	Devices and Printers	es and Printers	Boys
	C Lase of Access Center	File Lopic	🛔 Jack Mix Properties	× I-bit)	Player (32-bit)	
Levels	Fonts	lindesting	Current Luter Levels Advanced	s Settings	Graphics Settings	oner (Bealt) ck Andro
	finternet Options	🏄 Java (32	1 50 M			
FrontMic: 90	Mouse	Motwork Center	Minnahene Tard		r Options	(and a
Microphone Poacture 20 OdP	Programs and Postures	Recovery	a 0.02 • 10	/ Desktop	teApp and Desktop	
$\mathbf{M}(\mathbf{U}) = \mathbf{U}(\mathbf{U}) = U$	Security and Maintenance	Sound			ections be Seares	(500
	-				Providence .	300

7. If you still do not have a microphone signal, then Windows Microphone Privacy Setting may be Off.

B ....

Search for "Privacy Settings" and select Microphone. Switch on: Allow apps to access your microphone <On> Allow desktop apps to access your microphone <On>

Settings		– D ×
€ Home	Microphone	
Find a setting	Allow access to the microphone on this device	Know your privacy options
Privacy	If you allow access, people using this device will be able to choose if their apps have microphone access by using the settings on this page. Denying access blocks Windows features, Microsoft Store apps, and most desktop apps from accessing the microphone.	privacy Privacy dashboard Privacy dashboard
Windows permissions	Microphone access for this device is on	
B General	Change	Have a question?
위 Speech		Gernap
Inking & typing personalization	Allow apps to access your microphone If you allow access, you can choose which apps can access your	Make Windows better
Diagnostics & feedback	microphone by using the settings on this page. Denying access blocks apps from accessing your microphone.	Give us feedback
Activity history	On On	
pp permissions	Some desktop apps may still be able to access your microphone when settings on this page are off. Find out why	
5 Location	If an app is using your microphone, you'll see this icon: $\clubsuit$	
2 Camera		
] Microphone	Choose which Microsoft Store apps can access your microphone	
Voice activation	Some apps need to access your microphone to work as intended. Turning off an app here might limit what it can do.	
3 Notifications	3D Viewer On	
Account info	Camera On	
A Contacts		

8. You may need to close Scope and then open it again.